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
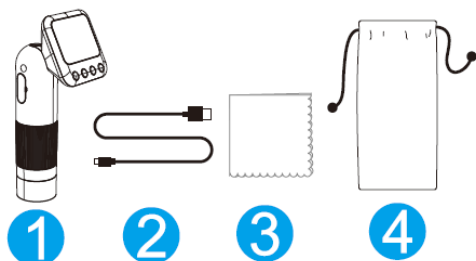
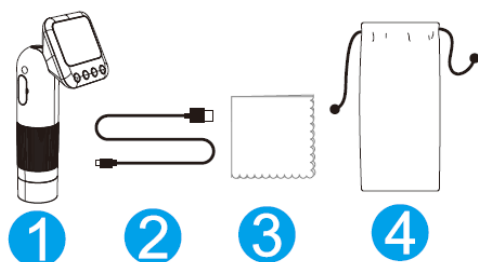
USER MANUAL – EN

IN APL-MS008CS Digital microscope Apexel 400X-800X with LCD display

SevenSport s.r.o. reserves the right to make any changes and improvements to its product without prior notice. Visit our website www.insportline.eu where you will find the latest version of the manual.

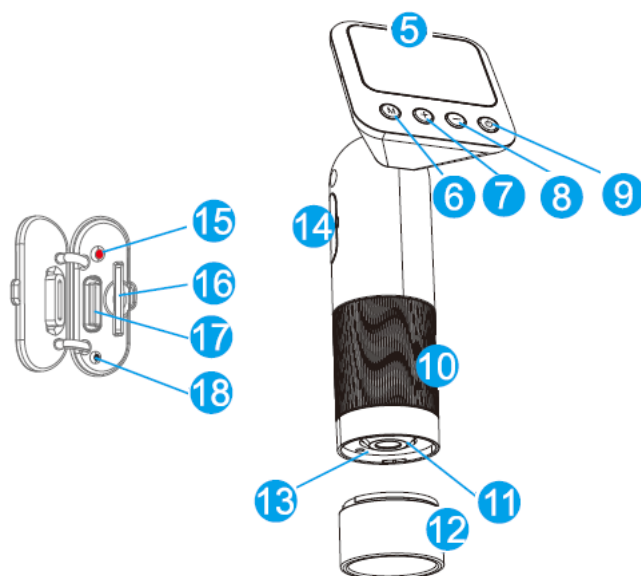1. Digital microscope
2. USB-C cable (data transfer, charging)
3. Cleaning cloth
4. Carrying bag

5. Display – display of the object (zoom 1x ~ 5x). You can save or record a screenshot of the display. Display of brightness and illumination.

6. M button

7. Button +

8. Button -

9. Power button

10. Focus – use the ring to focus on the subject

11. Lens – visually enlarges the object

12. Objective lens – used to track the object

13. LED ring – used to illuminate the monitored object

14. Output port – USB-C port for charging and data transfer. Cover the port with a cap when not in use.

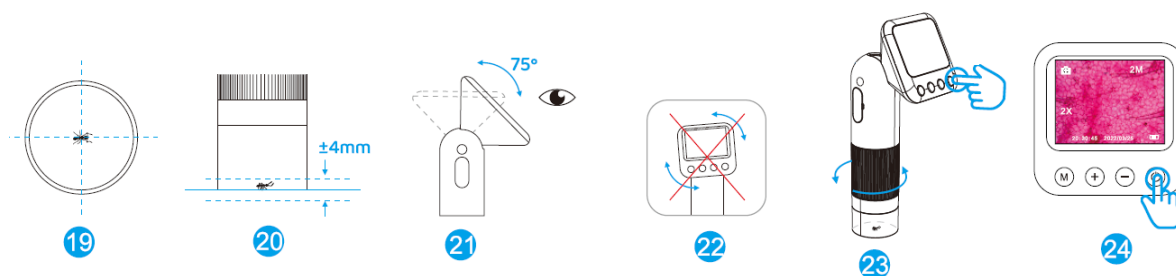
15. Charging indicator

16. TF card slot

17. USB-C port

18. Reset

USE

Microscope inspection

Hold down the button (9.) to start the display.

Hold down the button (7.) to start the light. Short press (7. / 8.) to change the zoom. Hold down the (8.) button to turn off the light.

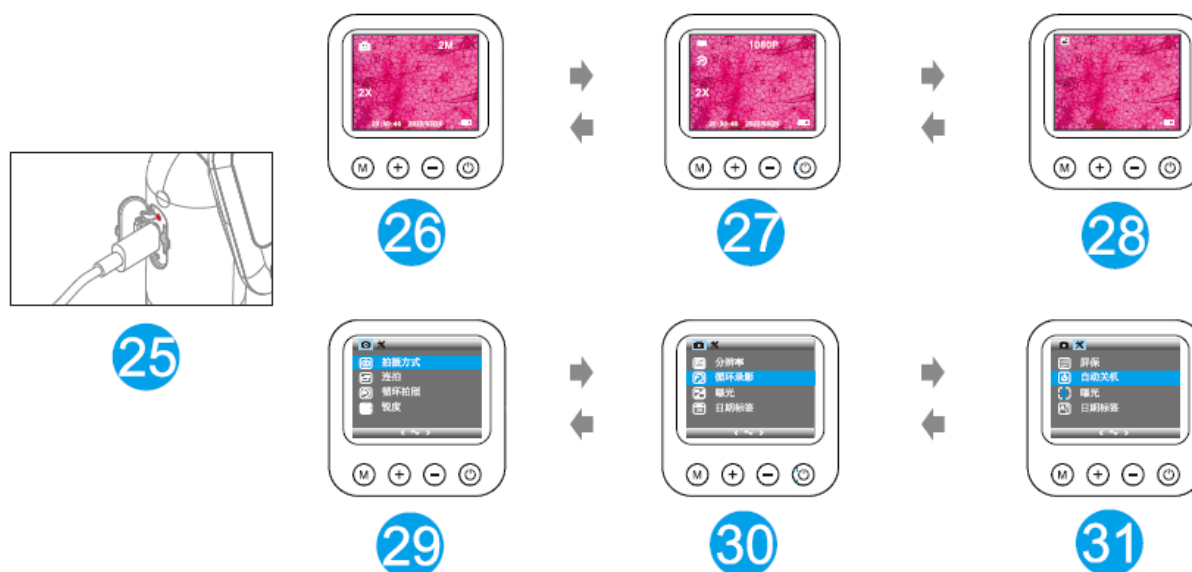
Start observing

Point the microscope at the object and make sure it is centered (19.). The height from the objective lens should be $\pm 4\text{mm}$ (20.).

Adjust the display angle to your needs (21.). The display can be positioned at an angle of 75° , otherwise you may damage the holder (22.).

Carefully focus the image and observe the changes. You can record a photo or video using the buttons (23. / 24.).

Controls

Press the power button (9.) for 2 seconds to start/stop the microscope.

Taking a photo/video and playing it back

Press the (6.) button to switch between modes (26. / 27. / 28.).

While taking a photo/video, press the button to zoom in (7.) or zoom out (8.), press the power button to take a photo/video.

Hold down (7.) to increase brightness, hold (8.) to decrease brightness. Hold the button (6.) to enter the setting (29. / 30. / 31.).

Playback of photos/videos

Press (7.) and (8.) to select pictures, or (9.) to play. During playback, press (7.) and (8.) to control the video.

Hold down the (6.) button to delete the file.

After entering the menu, press (6.) to switch between settings (photo / video / general), press (7.) and (8.) to cycle through the menu. To select the menu, press the power button. Press button (6.) to go back.

26. Photo interface

27. Video Interface

28. Playback

29. Photo settings interface

30. Video settings interface

31. Setup Interface

ADDITIONAL INFORMATION

Charging

Charge using the USB-C port and adapter. Charging time is approx. 2 hours, use time on a single charge is approx. 2 hours. While charging, the red LED lights up, after charging, the green LED lights up.

FAQ

Error	Cause	Solution
The microscope does not start	Low battery	Charge
The picture is not clear and is dark	The light is not on	Turn on the light
System stuck	Press the Restart button	Using a narrow object, press the restart button
The image is not sharp	Poorly focused, dirty surface	Check focus, clean
The display is too bright	Incorrectly set	Adjust the brightness in the settings

- Clean with soft and dry cloth.
- To connect to a computer, visit: www.apexellens.com
- Store in a shady and cool place.

ENVIRONMENT PROTECTION

After the product lifespan expires or if the possible repairing is uneconomical, dispose of it according to the local laws and is environmentally friendly in the nearest scrapyards.

By proper disposal you will protect the environment and natural resources. Moreover, you can help protect human health. If you are not sure of the correct disposing, ask local authorities to avoid law violations or sanctions.

Don't put the batteries among house waste but hand them in to the recycling place.

TERMS AND CONDITIONS OF WARRANTY, WARRANTY CLAIMS

General Conditions of Warranty and Definition of Terms

All Warranty Conditions stated here under determine Warranty Coverage and Warranty Claim Procedure. Conditions of Warranty and Warranty Claims are governed by Act No. 89/2012 Coll. Civil Code, and Act No. 634/1992 Coll., Consumer Protection, as amended, also in cases that are not specified by these Warranty rules.

The seller is SEVEN SPORT s.r.o. with its registered office in Kutnohorská 531, Štěrboholy Retail Park, 109 00, Prague 10, ČR, Company Registration Number: 26847264, registered in the Trade Register at Regional Court in Prague, Section C, Insert No. 116888.

According to valid legal regulations, it depends whether the Buyer is the End Customer or not.

"The Buyer who is the End Customer" or simply the "End Customer" is the legal entity that does not conclude and execute the Contract in order to run or promote his own trade or business activities.

"The Buyer who is not the End Customer" is a Businessman that buys Goods or uses services for the purpose of using the Goods or services for his own business activities. The Buyer conforms to the General Purchase Agreement and business conditions.

These Conditions of Warranty and Warranty Claims are an integral part of every Purchase Agreement made between the Seller and the Buyer. All Warranty Conditions are valid and binding, unless otherwise specified in the Purchase Agreement, in the Amendment to this Contract or in another written agreement.

Warranty Conditions

Warranty Period

The Seller provides the Buyer a 24 months Warranty for Goods Quality, unless otherwise specified in the Certificate of Warranty, Invoice, Bill of Delivery or other documents related to the Goods. The legal warranty period provided to the Consumer is not affected.

By the Warranty for Goods Quality, the Seller guarantees that the delivered Goods shall be, for a certain period of time, suitable for regular or contracted use, and that the Goods shall maintain its regular or contracted features.

Batteries

6-month battery warranty – we guarantee that battery's nominal capacity does not fall below 70% of its total capacity within 6 months of the product's sale.

The Warranty does not cover defects resulting from (if applicable):

- User's fault, i.e. product damage caused by unqualified repair work, improper assembly, insufficient insertion of seat post into frame, insufficient tightening of pedals and cranks
- Improper maintenance
- Mechanical damage
- Regular use (e.g. wearing out of rubber and plastic parts, moving mechanisms, joints, wear of brake pads/blocks, chain, tires, cassette/multi wheel etc.)
- Unavoidable event, natural disaster
- Adjustments made by unqualified person
- Improper maintenance, improper placement, damages caused by low or high temperature, water, inappropriate pressure, shocks, intentional changes in design or construction etc.

Warranty Claim Procedure

The Buyer is obliged to check the Goods delivered by the Seller immediately after taking the responsibility for the Goods and its damages, i.e. immediately after its delivery. The Buyer must check the Goods so that he discovers all the defects that can be discovered by such check.

When making a Warranty Claim the Buyer is obliged, on request of the Seller, to prove the purchase and validity of the claim by the Invoice or Bill of Delivery that includes the product's serial number, or eventually by the documents without the serial number. If the Buyer does not prove the validity of the Warranty Claim by these documents, the Seller has the right to reject the Warranty Claim.

If the Buyer gives notice of a defect that is not covered by the Warranty (e.g. in the case that the Warranty Conditions were not fulfilled or in the case of reporting the defect by mistake etc.), the Seller is eligible to require a compensation for all the costs arising from the repair. The cost shall be calculated according to the valid price list of services and transport costs.

If the Seller finds out (by testing) that the product is not damaged, the Warranty Claim is not accepted. The Seller reserves the right to claim a compensation for costs arising from the false Warranty Claim.

In case the Buyer makes a claim about the Goods that is legally covered by the Warranty provided by the Seller, the Seller shall fix the reported defects by means of repair or by the exchange of the damaged part or product for a new one. Based on the agreement of the Buyer, the Seller has the right

to exchange the defected Goods for a fully compatible Goods of the same or better technical characteristics. The Seller is entitled to choose the form of the Warranty Claim Procedures described in this paragraph.

The Seller shall settle the Warranty Claim within 30 days after the delivery of the defective Goods, unless a longer period has been agreed upon. The day when the repaired or exchanged Goods is handed over to the Buyer is considered to be the day of the Warranty Claim settlement. When the Seller is not able to settle the Warranty Claim within the agreed period due to the specific nature of the Goods defect, he and the Buyer shall make an agreement about an alternative solution. In case such agreement is not made, the Seller is obliged to provide the Buyer with a financial compensation in the form of a refund.

CZSEVEN SPORT s.r.o.

Registered Office: Kutnohorská 531, Štěrboholy Retail Park,
109 00, Prague 10, Czech Republic

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