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
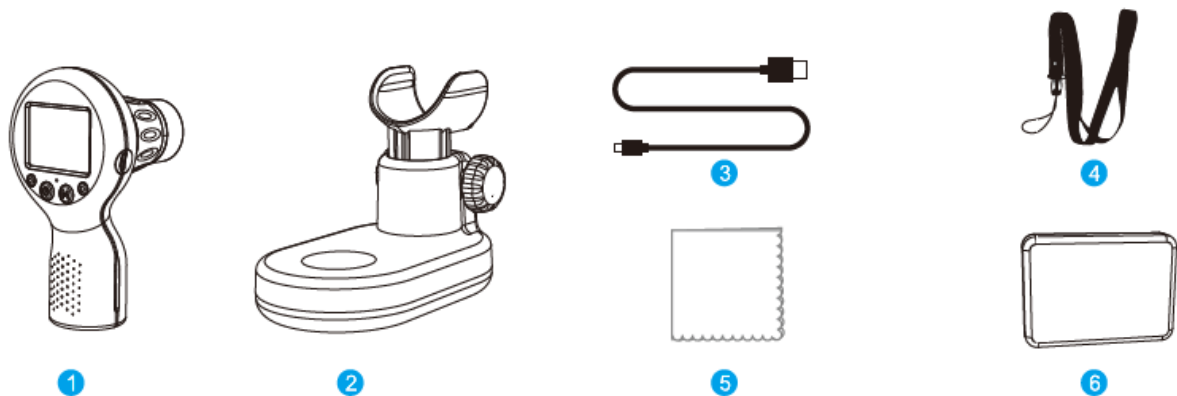
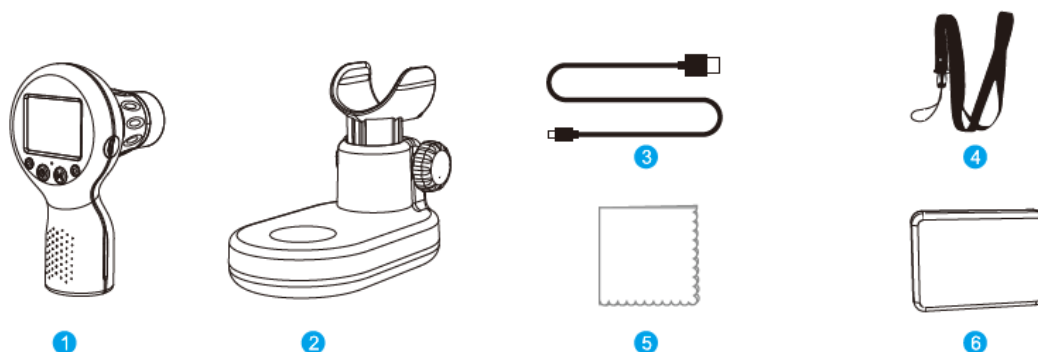
USER MANUAL – EN

IN APL-MS201 Children's digital microscope Apexel





SevenSport s.r.o. reserves the right to make any changes and improvements to its product without prior notice. Visit our website www.insportline.eu where you will find the latest version of the manual.

- | | |
|------------------------|-------------------|
| 1. Handheld microscope | 4. Hanging strap |
| 2. Microscope base | 5. Cleaning cloth |
| 3. USB-C cable | 6. Utensil box |

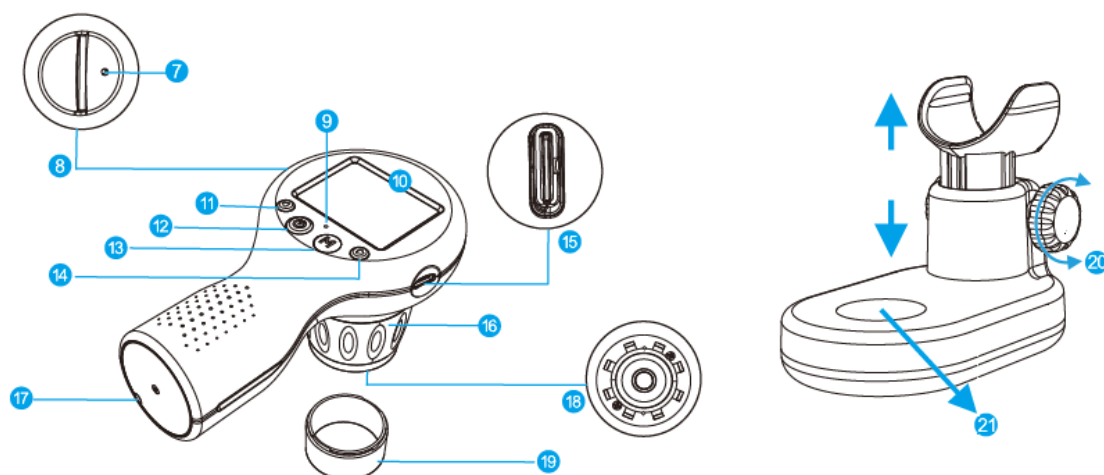7. Reset button	15. USB-C charging port
8. TF card slot	16. Focus ring
9. Charging indicator: red (low battery) / green (charged battery)	17. Strap hole
10. Display: 1x – 5x magnification viewing display. View object, display settings, take photo/video of viewed object	18. LED light – illumination of the monitored object
11. Decrease brightness (hold) / move up, digital zoom out (press)	19. Transparent cover
12. Power button (hold) / confirm, take photo (press)	20. Position adjustment knob
13. Menu button (hold) / switch between photos, record video, play (press)	21. Viewing area
14. Brightness Up (Hold) / Scroll Down, Digital zoom (Press)	

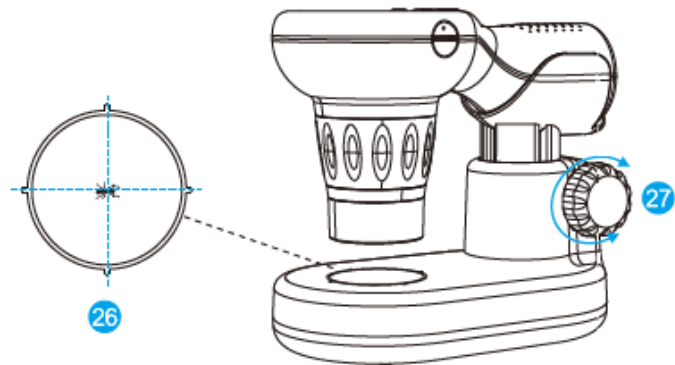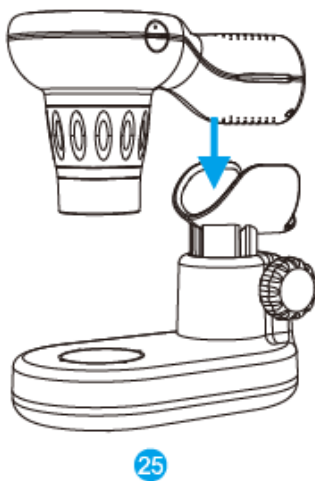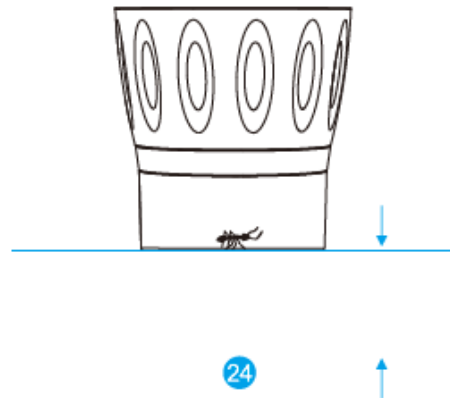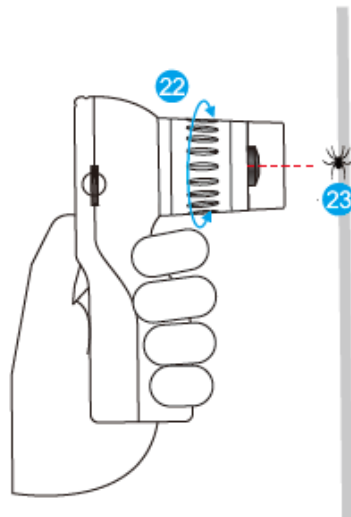1-1. Hold down the button (12) to start the microscope, the display will light up.

1-2. Press the button (11) to zoom out the image, press the button (14) to zoom in the image.

2-1. Place the tracked object in the center of the lens (23) and use the focus ring (22) to focus. Use the buttons (11 and 14) to digitally zoom in or zoom out the image. Press the button (12) to take a photo or record a video.

3-1. You can place the microscope in the holder (25).

3-2. Place the tracked object in the center of the viewing area (26).

3-3. Use the knob (27) to adjust the height of the microscope base and use the ring (22) to adjust the focus.

Notes:

- Make sure the microscope battery is sufficiently charged before use.
- When tracking an object, we recommend focusing using the focus ring and then using the zoom function.
- Maintain a suitable distance and viewing angle.
- Do not move the microscope while taking a photo or video.
- Always turn off the microscope after use and clean the lens and viewing area with a cloth.

1. Switch on: hold down the button (12) for 2 seconds to start.
2. Switch off: hold the button (12) for 2 seconds to switch off.
3. To take a photo / video / change mode: press the button (13) to switch between the interface (28 / 29 / 30).
4. Taking a photo/video: while taking a photo or video, you can use the buttons (11 and 14) to digitally zoom in or zoom out the image. Press the button (12) to take a photo or video.
5. To start the light / adjust the brightness: hold the button (14) to start the light or increase the brightness, hold the button (11) to decrease the brightness or turn off the light.
6. Setting: hold the button (13) to enter the setting interface. Use the buttons (11 / 14) to navigate the settings. Press the button (12) to enter the selected setting.
7. Adjust settings: use the buttons (11 / 14) to adjust the selected value. Press button (12) to confirm the setting, press button (13) to return to the setting.
8. To view photos/videos: press (11 / 14) to browse the files, press (12) to select a file.
9. Video playback: hold down the buttons (11/14) to fast-forward or rewind the video.
10. To delete a file: to delete a file, hold down the button (13).

28. Photo interface

29. Video Interface

30. Display interface

31. Delete a video

ADDITIONAL INFORMATION

Charging

Charge using the USB-C port and adapter. Charging time is approx. 1.5 hours, use time on a single charge is approx. 3 hours. While charging, the red LED lights up, after charging, the green LED lights up.

Insert memory card back side facing up.

FAQ

Error	Cause	Solution
The microscope does not start	Low battery	Charge
The picture is not clear and is dark	The light is not on	Turn on the light
System stuck	Press the Restart button	Using a narrow object, press the restart button
The image is not sharp	Poorly focused, dirty surface	Check focus, clean
The display is too bright	Incorrectly set	Adjust the brightness in the settings

Precautions

- Do not expose to direct sunlight.
- Protect lenses against dirt and damage.

- Do not place the microscope on an uneven surface.
- To protect your eyes, limit the length of use.

Notification

- Clean with a soft and dry cloth.
- To connect to a computer, visit: www.apexellens.com
- Store in a shady and cool place.

ENVIRONMENT PROTECTION

After the product lifespan expires or if the possible repairing is uneconomical, dispose of it according to the local laws and is environmentally friendly in the nearest scrapyard.

By proper disposal you will protect the environment and natural resources. Moreover, you can help protect human health. If you are not sure of the correct disposing, ask local authorities to avoid law violations or sanctions.

Don't put the batteries among house waste but hand them in to the recycling place.

TERMS AND CONDITIONS OF WARRANTY, WARRANTY CLAIMS

General Conditions of Warranty and Definition of Terms

All Warranty Conditions stated here under determine Warranty Coverage and Warranty Claim Procedure. Conditions of Warranty and Warranty Claims are governed by Act No. 89/2012 Coll. Civil Code, and Act No. 634/1992 Coll., Consumer Protection, as amended, also in cases that are not specified by these Warranty rules.

The seller is SEVEN SPORT s.r.o. with its registered office in Kutnohorská 531, Štěrboholy Retail Park, 109 00, Prague 10, ČR, Company Registration Number: 26847264, registered in the Trade Register at Regional Court in Prague, Section C, Insert No. 116888.

According to valid legal regulations, it depends whether the Buyer is the End Customer or not.

"The Buyer who is the End Customer" or simply the "End Customer" is the legal entity that does not conclude and execute the Contract in order to run or promote his own trade or business activities.

"The Buyer who is not the End Customer" is a Businessman that buys Goods or uses services for the purpose of using the Goods or services for his own business activities. The Buyer conforms to the General Purchase Agreement and business conditions.

These Conditions of Warranty and Warranty Claims are an integral part of every Purchase Agreement made between the Seller and the Buyer. All Warranty Conditions are valid and binding, unless otherwise specified in the Purchase Agreement, in the Amendment to this Contract or in another written agreement.

Warranty Conditions

Warranty Period

The Seller provides the Buyer a 24 months Warranty for Goods Quality, unless otherwise specified in the Certificate of Warranty, Invoice, Bill of Delivery or other documents related to the Goods. The legal warranty period provided to the Consumer is not affected.

By the Warranty for Goods Quality, the Seller guarantees that the delivered Goods shall be, for a certain period of time, suitable for regular or contracted use, and that the Goods shall maintain its regular or contracted features.

Batteries

6-month battery warranty – we guarantee that battery's nominal capacity does not fall below 70% of its total capacity within 6 months of the product's sale.

The Warranty does not cover defects resulting from (if applicable):

- User's fault, i.e. product damage caused by unqualified repair work, improper assembly, insufficient insertion of seat post into frame, insufficient tightening of pedals and cranks

- Improper maintenance
- Mechanical damage
- Regular use (e.g. wearing out of rubber and plastic parts, moving mechanisms, joints, wear of brake pads/blocks, chain, tires, cassette/multi wheel etc.)
- Unavoidable event, natural disaster
- Adjustments made by unqualified person
- Improper maintenance, improper placement, damages caused by low or high temperature, water, inappropriate pressure, shocks, intentional changes in design or construction etc.

Warranty Claim Procedure

The Buyer is obliged to check the Goods delivered by the Seller immediately after taking the responsibility for the Goods and its damages, i.e. immediately after its delivery. The Buyer must check the Goods so that he discovers all the defects that can be discovered by such check.

When making a Warranty Claim the Buyer is obliged, on request of the Seller, to prove the purchase and validity of the claim by the Invoice or Bill of Delivery that includes the product's serial number, or eventually by the documents without the serial number. If the Buyer does not prove the validity of the Warranty Claim by these documents, the Seller has the right to reject the Warranty Claim.

If the Buyer gives notice of a defect that is not covered by the Warranty (e.g. in the case that the Warranty Conditions were not fulfilled or in the case of reporting the defect by mistake etc.), the Seller is eligible to require a compensation for all the costs arising from the repair. The cost shall be calculated according to the valid price list of services and transport costs.

If the Seller finds out (by testing) that the product is not damaged, the Warranty Claim is not accepted. The Seller reserves the right to claim a compensation for costs arising from the false Warranty Claim.

In case the Buyer makes a claim about the Goods that is legally covered by the Warranty provided by the Seller, the Seller shall fix the reported defects by means of repair or by the exchange of the damaged part or product for a new one. Based on the agreement of the Buyer, the Seller has the right to exchange the defected Goods for a fully compatible Goods of the same or better technical characteristics. The Seller is entitled to choose the form of the Warranty Claim Procedures described in this paragraph.

The Seller shall settle the Warranty Claim within 30 days after the delivery of the defective Goods, unless a longer period has been agreed upon. The day when the repaired or exchanged Goods is handed over to the Buyer is considered to be the day of the Warranty Claim settlement. When the Seller is not able to settle the Warranty Claim within the agreed period due to the specific nature of the Goods defect, he and the Buyer shall make an agreement about an alternative solution. In case such agreement is not made, the Seller is obliged to provide the Buyer with a financial compensation in the form of a refund.

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